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

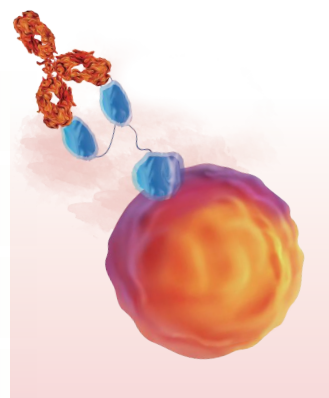
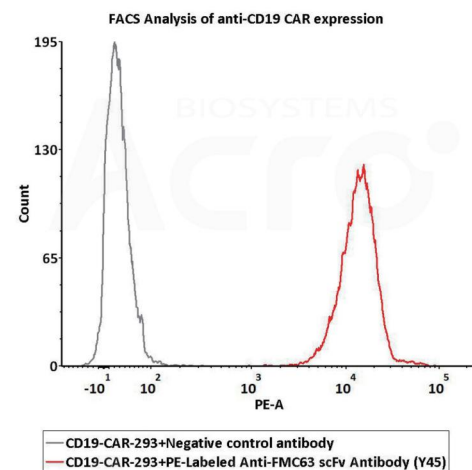
5e5 of the Anti-CD19 CAR-293 cells were stained with 100 µL of 1:50 dilution (2 µL stock solution in 100 µL FACS buffer) of PE-Labeled Monoclonal Anti-FMC63 scFv Antibody, Mouse IgG1 (Cat. No. FM3-HPY53) and negative control antibody (Cat. No. DNP-PM1) respectively, PE signal was used to evaluate the binding activity (QC tested).

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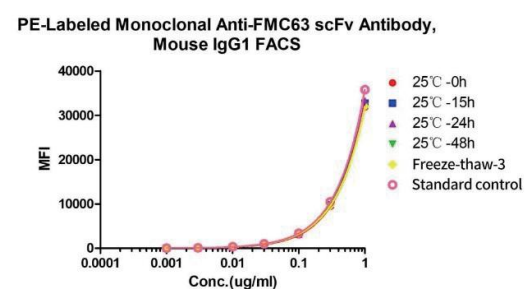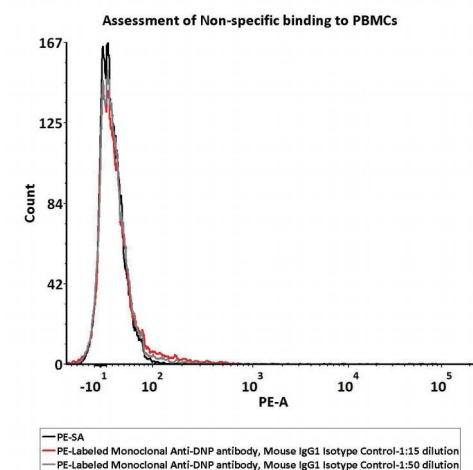
PE-Labeled Monoclonal Anti-FMC63 scFv Antibody, Mouse IgG1 (Cat. No. FM3-HPY53) at 25°C for 48 hours and freeze-throw cycles are stable without performance reduction.

5e5 of PBMCs were stained with a series of concentrations of PE-Labeled Monoclonal Mouse IgG1 Isotype Control (Cat. No. DNP-PM1), and PE signal was used to evaluate the binding activity (QC tested).

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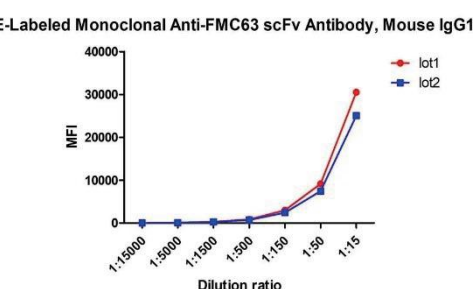Binding activity of two different lots of PE-Labeled Monoclonal Anti-FMC63 scFv Antibody, Mouse IgG1 (Cat. No. FM3-HPY53) against Anti-CD19 CAR-293 cells was evaluated by flow cytometry. The result shows very high batch-to-batch consistency.

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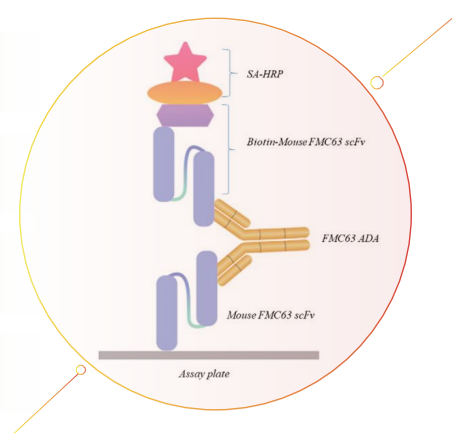
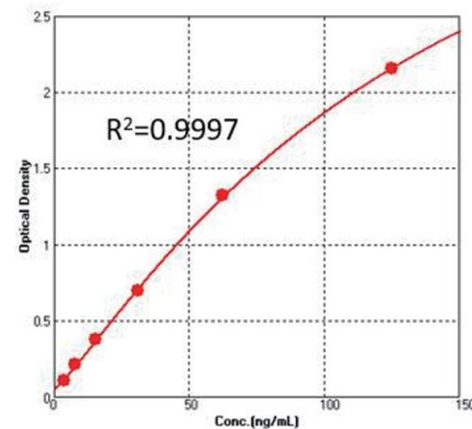
Detection of FMC63 ADA titer by bridging-ELISA Assay. Immobilized Mouse FMC63 scFv can bind FMC63 ADA in 1:10 human serum, and then add Biotin-Mouse FMC63 scFv. Detection was performed using HRP-conjugated streptavidin with sensitivity of 78 ng/mL (QC tested).

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アッセイの種類	製品番号	ホスト	説明
ADA Assay	FM3-S93	Rabbit	Rabbit Anti-Mouse FMC63 scFv Polyclonal Antibody (recommended for ADA assay) (MALS verified)
	CD9-M52Hb	Mouse	FMC63 scFv, His Tag (SPR verified)
	CD9-BV4324b	Mouse	Biotinylated Anti-CD19 scFv, Fc, Avitag™ (FMC63) (MALS verified)
Neutralizing Assay	FM3-Y45A1	Mouse	Monoclonal Anti-FMC63 scFv Antibody, Mouse IgG1 (Y45) (Carrier-free) (recommended for ADA assay) DMF Filed

Her2 BAFFR LAG-3  
Fc Receptor Siglec-10  
Biotinylated Protein  
PD-L1 VEGF165 CD3 epsilon  
CD20 PD-1 BCMA  
CD27 PVRIG  
CD47 PSMA  
CFGL1 TFPI  
Siglec-15 Integrin  
CD24 CD3E & CD3D CD20  
CD19 FcRn PCSK9  
IL-2 R alpha  
CAR-T Target Protein  
Glypican 3 Integrin  
ADA Service  
EGF R B7-H3 BCMA  
Integrin TIGIT TGF-beta 1  
CD200 GITR Nectin-4  
CD20 VEGF165 CD3D FGLI  
CD40 CD137 CD137L  
BCMA CD137L  
SIRP alpha ADA Service  
Nectin-4 Biotinylated Protein CD3E & CD3D  
SPR /BLI analytical service

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